

“Special K” Trace Elements Recipe devised by Merchant group (Janette Kropat and Davin Malasarn) – work is unpublished, details of development and analysis will follow. Please do not distribute outside the NAABB consortium until after publication.

Revised June 2, 2010

Make preliminary concentrated stock solutions in Part A first, and, where indicated, use these to make the individual stock solutions in Part B listed below. Only solutions in Part B are added directly to media.

A. Preliminary concentrated stock solutions

Pre-1. EDTA-Na ₂ concentrate	125 mM	13.959 g in ~ 250 ml, titrate to pH 8.0 with trace element grade KOH (~1.7 g), and bring up to a volume of 300 ml
Pre-2. (NH ₄) ₆ Mo ₇ O ₂₄ concentrate	285 μM	(NH ₄) ₆ Mo ₇ O ₂₄ ·4H ₂ O: 0.088 g, bring up to a volume of 250 mL
Pre-3. Na ₂ SeO ₃ concentrate	1 mM	Na ₂ SeO ₃ : 0.043 g, bring up to a volume of 250 mL

B. Individual Stock Solutions for medium (1000×)

Bring each stock solution up to 250 mL in water. Use 1 mL of each individual stock solution in 1 L medium.

Stock Solution	Concentration in stock	Composition
1. EDTA-Na ₂	25 mM	EDTA-Na ₂ : 50 mL of 125 mM EDTA-Na ₂ concentrate (Pre-1) from Step A
2. (NH ₄) ₆ Mo ₇ O ₂₄	28.5 μM*	(NH ₄) ₆ Mo ₇ O ₂₄ ·4H ₂ O: 25 mL of 285 μM (NH ₄) ₆ Mo ₇ O ₂₄ concentrate (Pre-2) from Step A
3. Na ₂ SeO ₃	0.1 mM	Na ₂ SeO ₃ : 25 mL of 1 mM Na ₂ SeO ₃ concentrate (Pre-3) from Step A
4. Zn-EDTA	2.5 mM 2.75 mM	ZnSO ₄ ·7H ₂ O: 0.18 g EDTA-Na ₂ : 5.5 mL of 125 mM EDTA-Na ₂ concentrate (Pre-1) from Step A
5. Mn-EDTA	6 mM 6 mM	MnCl ₂ ·4H ₂ O: 0.297 g EDTA-Na ₂ : 12 mL of 125 mM EDTA-Na ₂ concentrate (Pre-1) from Step A

6. Fe·EDTA 20 mM FeCl₃·6H₂O: 1.35 g
 22 mM EDTA-Na₂: 2.05 g
 22 mM Na₂CO₃ (sodium carbonate): 0.58 g
(Combine EDTA-Na₂ with sodium carbonate in water and mix. Add FeCl₃·6H₂O after the first two components dissolve. Do Not Use Pre-1.)
7. Cu·EDTA 2 mM CuCl₂·2H₂O: 0.085 g
 2 mM EDTA-Na₂: 4 mL of 125 mM EDTA-Na₂ concentrate
(Pre-1) from Step A

Notes:

*The final [Mo] in the 1× medium is 0.2 μM

Total [EDTA] in 1× medium: (25 + 2.75 + 6 + 22 + 2) = 57.75 μM

Merchant lab
unpublished